



MycoFog® Quick Start Guide

Please consult full manual at <https://mycofog.com>

Day before use: Charge MycoFog® Device with supplied cable.



> 8 hours

Day of use: Take out all cell cultures from incubator, set Incubator to **37°C** or **45°C**, turn off gas control and remove water pan or turn off active humidification.

1. Open MycoFog®



Twist then
lift top



2. Pour MycoFog® Reagent into reagent reservoir of MycoFog®



3. Remove the wick holder, insert wick, replace holder and press firmly. Replace top of MycoFog®.




4. Place MycoFog® in incubator, mid-chamber, wait 5 minutes.



5 minutes

5. Press Start

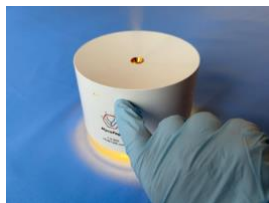

6. Close incubator door set timer.



180 Minutes (or **240** mins if using 37 °C)

7. Remove MycoFog®, if necessary return temperature to 37°C. Your Incubator is now decontaminated and ready for use!